


GUIDE TO DONNING AND DOFFING OF NON-STERILE CLEANROOM GLOVES

DONNING GLOVES

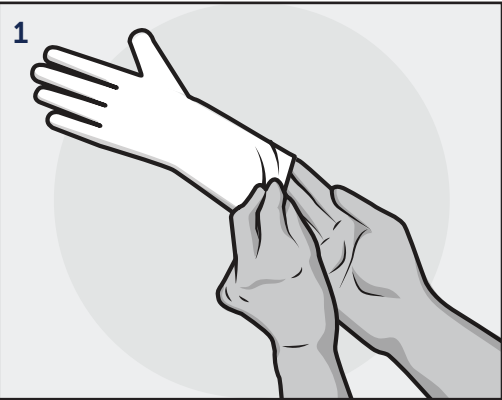


Grasp the beaded cuff of one glove and carefully pull over the opposite hand.




Gently ease the glove up over the hand and wrist.



Grasp the second glove at the edge of the beaded cuff.

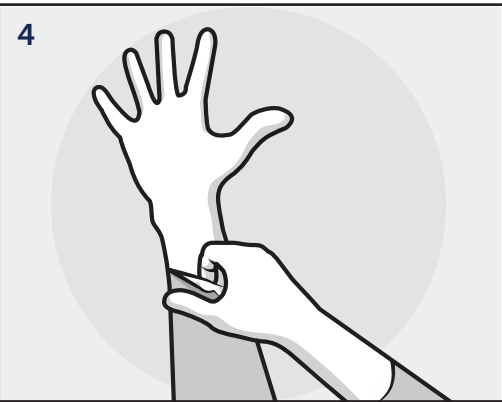


Gently ease the glove up over the hand and wrist.

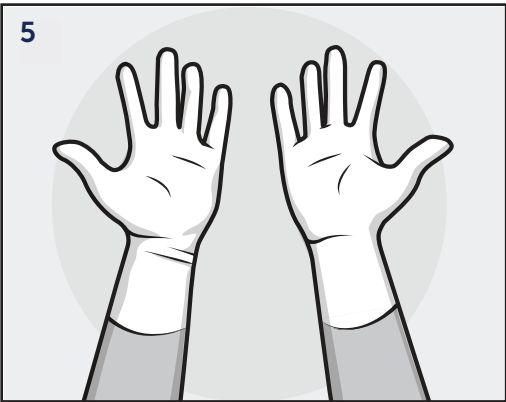


You are now ready to perform your task.


DOFFING GLOVES

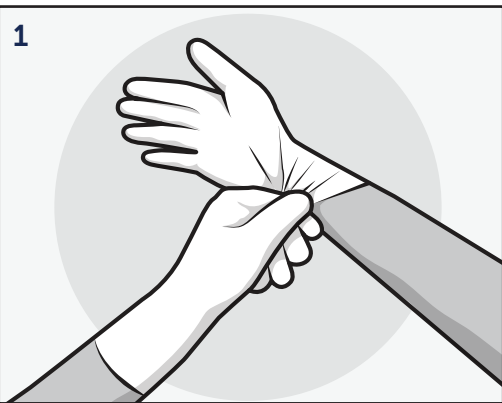

Grasp the exterior of one glove with your other gloved hand.

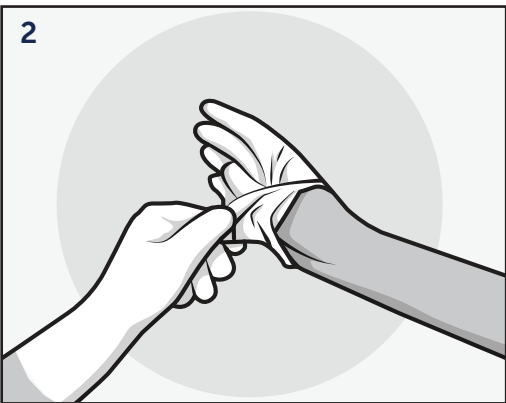


Carefully pull the glove off your hand, turning it inside-out.

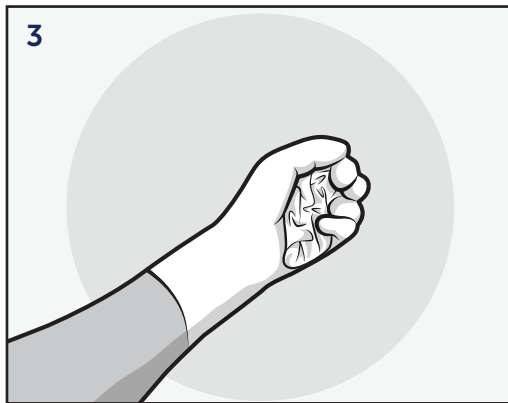


Ball the glove up and hold in your other gloved hand.



Slide your ungloved fingers into the opening of the other glove.

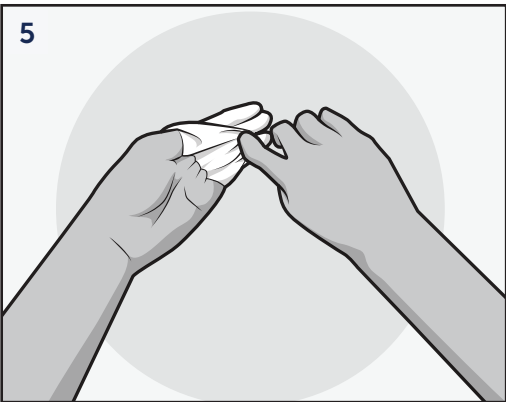

Carefully pull the glove off your hand, turning it inside-out and wrapping it around the doffed glove in your palm.

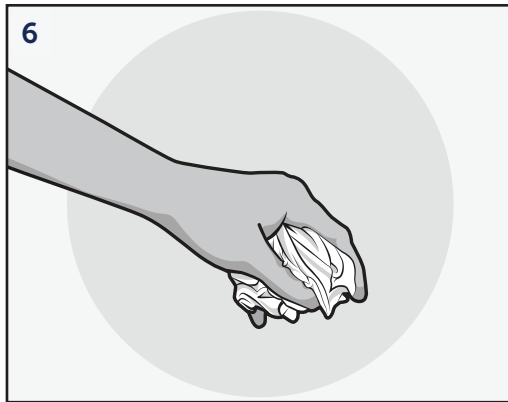

Discard appropriately.

For more information, contact your HALYARD* representative or distributor, or visit www.purezerogloves.com