



GUIDE TO DONNING AND DOFFING OF STERILE CLEANROOM GLOVES


DONNING GLOVES

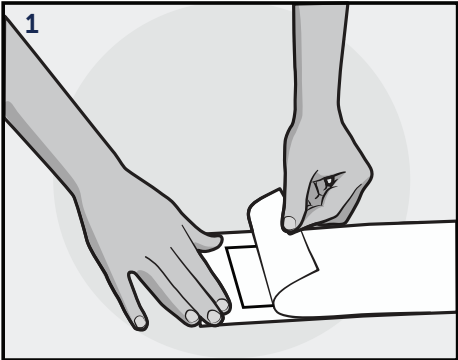


Peel open the pouch which keeps the product sterile. Open clear wallet.

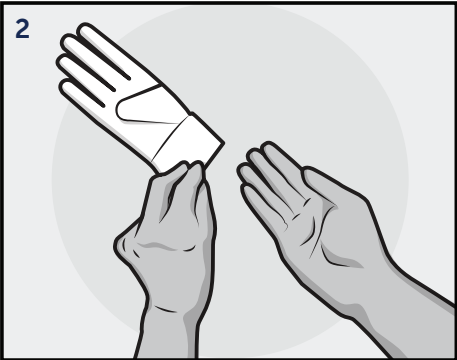

Grasp the interior of one glove and carefully pull over the opposite hand.

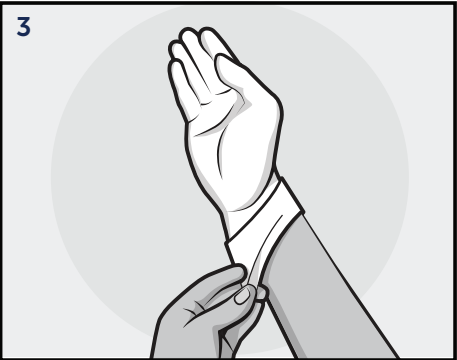


Gently ease the glove up over the hand and wrist, leaving the cuff folded.

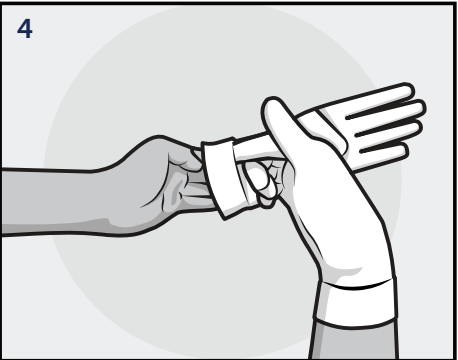


Lift the second glove under the folded cuff using gloved fingers.



Slide your second hand into the glove by holding it under the folded cuff, then unfold the cuff over the wrist and arm.

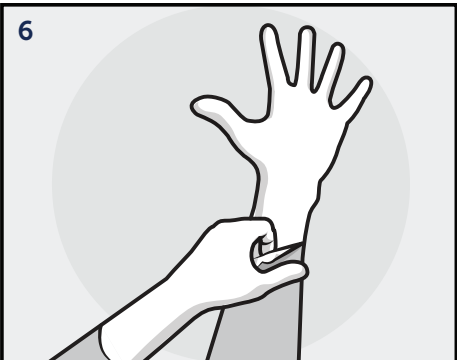


Using gloved fingers, unfold the second cuff over your wrist and arm.

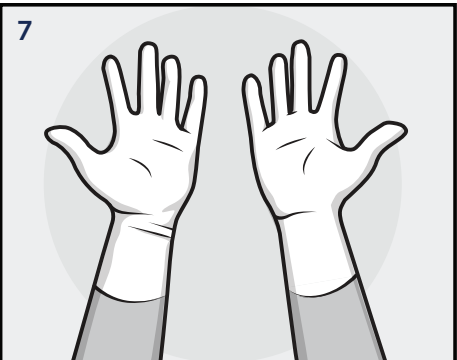


To remain sterile, do not touch anything with your gloved hands.


DOFFING GLOVES

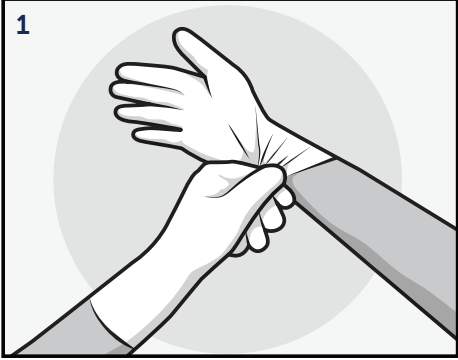


Grasp the exterior of one glove with your other gloved hand.

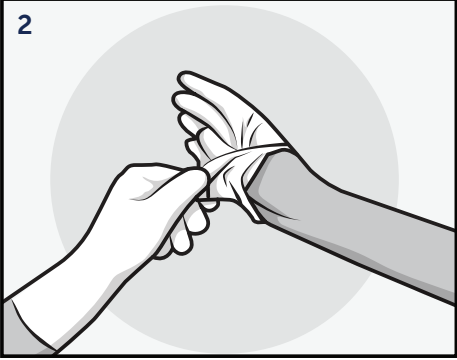


Carefully pull the glove off your hand, turning it inside-out.

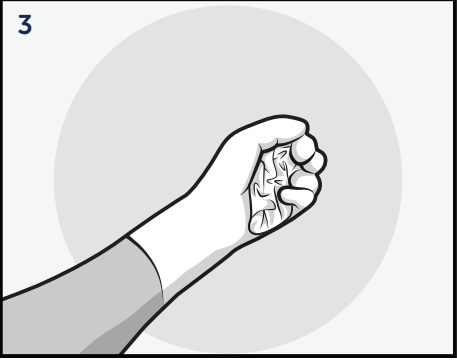


Ball the glove up and hold in your other gloved hand.

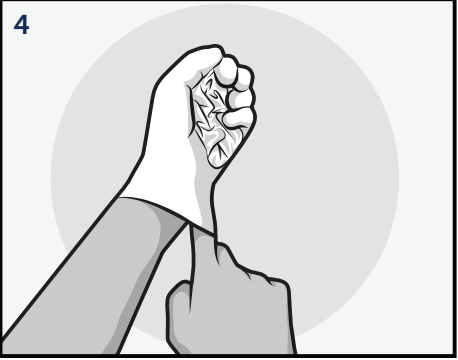

Slide your ungloved fingers into the opening of the other glove.

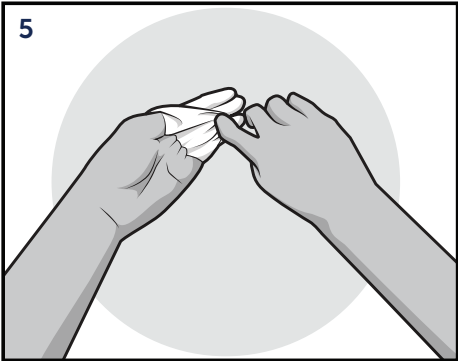

Carefully pull the glove off your hand, turning it inside-out and wrapping it around the doffed glove in your palm.

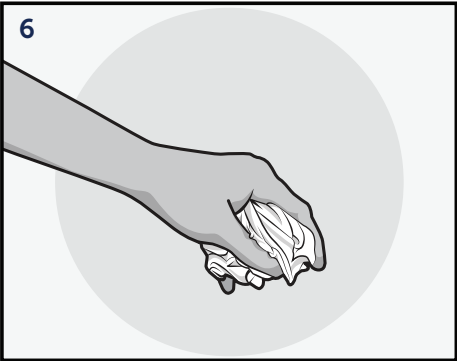

Discard appropriately.

HALYARD* recommends double donning/gloving. Don the gloves so that the sleeve cuffs are between the two pairs of gloves.

For more information, contact your HALYARD* representative or distributor, or visit www.purezerogloves.com